

UV-3200

Double Beam Spectrophotometer



LABINDIA®
Where commitment is culture





UV-3200 Double Beam Spectrophotometer

The two detectors measure sample and reference respectively and simultaneously for optimizing measurement accuracy. They provide excellent performance for measurements in the range of 190 to 1100nm. They are suitable for pharmaceutical, biochemical and clinical lab applications as well as routine applications such as quantitative analysis, kinetics, wavelength scan, multiple components and

DNA/Protein, PC Windows application software make these instruments versatile.

variables like (bandwidths) For Stand-alone instrument, all software methods are included as built-in standard; this eliminates the need of software. Online software upgrade via internet helps to keep it updated. Data Download-to-PC software expands the data storage to unlimited.

Specifications	Double Beam Optical System Model 3200
Wavelength Range	190nm to 1100nm
Working Mode	Stand alone and PC controlled with window based application software UV/VIS Analyst
Spectral Band Width	Variable 0.5,1,2,4 nm
Monochromator	Double beam hollow graphic gratings 1200 lines/mm
Wavelength Display	0.1nm
Wavelength setting	0.1nm
Wavelength accuracy	+ 0.1 nm @656.1nm D2 +0.3nm((190 to 1100nm)
Wavelength Repeatability	+0.1nm
Stray Light	0.02% @ 340nm for NaNO ₂ 0.25% @ 198nm for KCl
Photometric Range (Approx.)	0-200%T Absorbance -0.4 to 4
Photometric Accuracy	0.002 Abs(0.5) 0.002 Abs(1.0) 0.002 Abs(2.0)
Photometric Reproducibility	0.001 Abs(0.5 Abs) 0.001 Abs(1.0 Abs) 0.001 Abs(2.0 Abs)
Baseline Stability	< 0.0002 Abs/H @ 500 nm
Baseline Flatness	+ 0.0005 Abs
Noise Level	0.000016 Abs RMS @ 500 nm
Standard Cell Holder	Standard 10mm Single Cell Holder 2pcs
Light Source	Tungsten and Deuterium Lamp, pre-aligned
Output	USB Port and Parallel Port (Printer)
DNA/Protein Measurement	Included
Dimension (W x D x H)	600 x 450 x 200 mm
Weight	22kg
* Optional 8-cell changer	

The PC Models come standard with windows based application software UV/Vis Analyst


V-1100D/UV-1100D UV-3200 Local Control Software

All methods are included as built-in standard; this eliminates the need of software. Online software update via Internet . The local control software includes functions such as: Photometry, Quantitative, Wavelength Scan, Kinetics, DNA/Protein, Multi-wavelength Test and System Utilities.



Standard Curve


Up to 10 standard solutions may be used to establish calibration equation curve. There is a choice of four methods for fitting curve through the calibration points: Linear fit, Linear fit through zero, square fit and cubic fit.

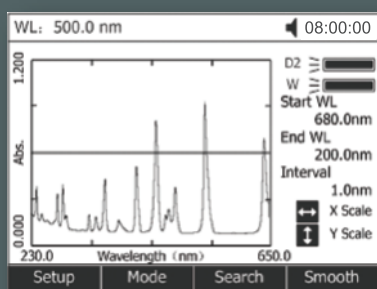

Wavelength Scan


The Wavelength Scan intervals are 0.1, 0.2, 0.5, 1, 2, 5nm, and High, Medium and Low scan speeds are available. Scan speeds vary from 100 to 1000 nm/min. Wavelengths are scanned from high to low so that the instrument stand-by at high wavelength. This minimizes the degradation of UV sensitive samples. Precise control of filter and lamp changes means that their effects are not seen on the final scan. Post-run manipulation includes re-scaling axes, curve tracking and peak picking.

V-1100D/UV-1100D

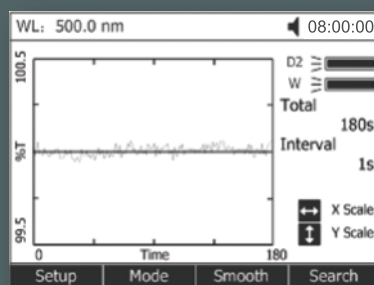


Kinetics

This mode may be used for time course scanning or reaction rate calculations. Abs. vs. time graphs is displayed on the screen in real time. Wait time and measurement time up to 12 hours may be entered with time intervals of 0.5, 1, 2, 5, 10, 30, seconds and 1 min. Post-run manipulation includes re-scaling, curve tracking and selection of the part of the curve required for the rate calculation. Rate is calculated using a linear regression algorithm before multiplying by the entered factor.

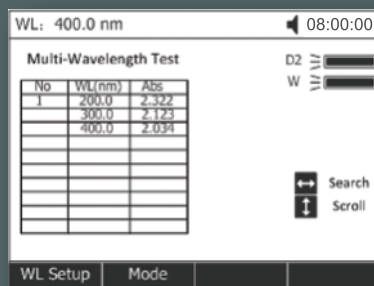

Multi-Wavelength

Up to 10 wavelengths may be entered, allowing the measurement of multiple wavelengths on a series of Samples.

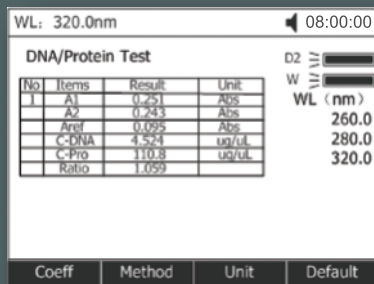

DNA/Protein Test

Concentration and DNA purity are calculated
Absorbance ratios 260nm/280nm or 260nm/230nm with optional subtracted absorbance at 320nm
DNA Concentration= $62.9 \cdot A_{260} - 36.0 \cdot A_{280}$
Or $49.1 \cdot A_{260} - 3.48 \cdot A_{230}$
Protein Concentration= $1552 \cdot A_{260} - 757.3 \cdot A_{280}$
Or $183 \cdot A_{260} - 75.8 \cdot A_{230}$
Other wavelengths and factors may be entered.


PC Application software

The Labindia Windows®

based PC application software UV/Vis Analyst takes the best features of the stand-alone version plus more powerful data processing, expanded data collecting, and storage capability. It comes standard with MAPADA PC models and is optional to stand-alone models.



The PC application software offers:

1. Photometric Mode
2. Quantitative test(standard curve)
3. Wavelength Scan
4. Kinetics
5. DNA/Protein
6. Multi-Wavelength
7. System Utility

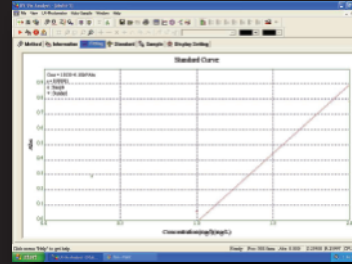

● Kinetics (Abs vs. Time)

The Kinetics mode may be used for time course scanning or reaction rate calculations. Abs.Vs.Time graphs are displayed on the screen in real time. Waiting time, measurement time and time intervals may be entered. Post-run manipulation includes re scaling, curve tracking and selection of the part of the curve required for the rate calculation. Rate is calculated using a linear regression algorithm before multiplying by the entered factor.

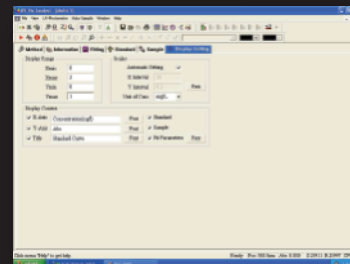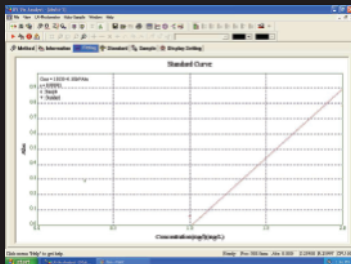

● Quantitative Test (Standard curve)

Use up to 20 standards to establish standard curve. Four methods for fitting a curve:

1. Linear fit
2. Linear through zero
3. Square fit
4. Cubic fit



● DNA/Protein

Concentration and DNA purity are quickly and easily calculated: Absorbance ratios 260nm/280nm with optional subtracted absorbance at 320nm.

DNA Concentration= $62.9 \times A_{260} - 36.0 \times A_{280}$

Protein Concentration= $1552 \times A_{260} - 757.3 \times A_{280}$

Other wavelengths and factors may be entered.

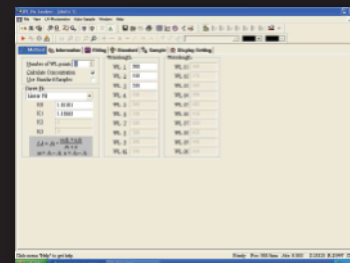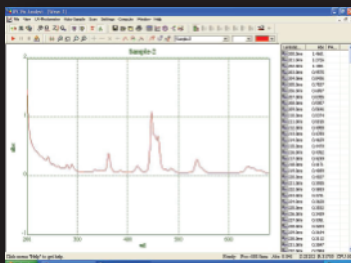

● Wavelength Scan

Automatically record peaks and valleys. The quantity of channels is unlimited; you can simultaneously store as many as desired. Post-run manipulation and processing includes:

1. Re-scaling axes, curve
2. 1st to 4th derivative
3. Smoothing, combination, zooming, overlap.


● Multi-wavelength

Up to 20 wavelengths can be selected and multiple samples can be measured. (Auto cell changer is required to run multiple samples automatically)

Optional Accessories


Micro Cell Holder

900210

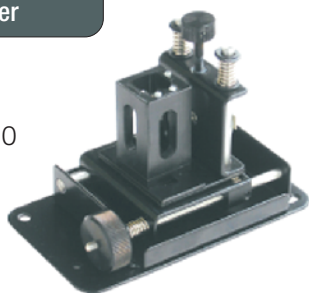

Square Cuvettes Glass

Square Cuvettes Quarts
916111-916115

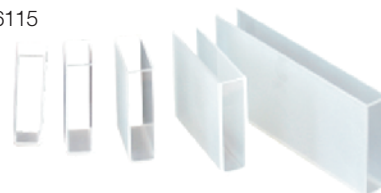

Test Tube Holder

900530



Mirco Cell Quartz

916121-916123

Self Masking Cont. Flowthrough G.Cell

916131-916134

Self Masking Cont. Flowthrough Q.Cell

916141-916144



8-Position Auto Cell Changer

900310



Stylus Printer

900920



Solid Sample Holder (Single Cell)

900550




Milas Deuterium Lamp

916633



Water-Jacketed Cell Holder

900610

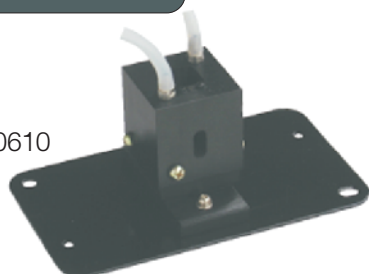

Constant-Temperature Sipper System

900130

Constant-Temperature System
900120

Sipper System
900110



LABINDIA®

Where commitment is culture

www.labindia.com / For Enquiries: sales.mfd@labindia.com

Labindia Analytical Instruments Pvt. Ltd.

FACTORY: Plot No. EL-72, Electronic Zone, TTC Industrial Area, Thane Belapur Road, Navi Mumbai - 400 705.
Tel.: +91-22-2762 6660/61/62/63/65163480/81
Fax: +91-22-27626664 E-mail: lipfactory@vsnl.net

APPLICATION & TRAINING CENTRE:

R-909, TTC Industrial Area, Thane Belapur Road, Rabale, Navi Mumbai - 400 701. Tel.: +91-22-2760 6955
Fax: +91-22-2760 6706

HEAD OFFICE & REGIONAL OFFICES:

Thane: 201, Nand Chambers, LBS Marg, Near Vandana Cinema, Thane (W) - 400 602. Tel.: +91-22-2598 6075 / 2598 6000
Fax: +91-22-2541 0420 / 2533 5940 E-mail: sales.mfd@labindia.com
Delhi: G-4, Pal Mohan Sadan, 26/32, East Patel Nagar, New Delhi - 110 008. Tel.: +91-11-4330 6001 / 10
Fax: +91-11-2585 1066
Chennai: B-1, Alsa Regency, 165, Eldams Road, Chennai - 600 018.
Tel.: +91-44-2434 7008 / 2432 0352 Fax: +91-44-2434 6328
Kolkatta: 165-A, S. P. Mukherjee Road, Kolkatta - 700 026.
Tel.: +91-33-2466 1396 / 2466 3362 Fax: +91-11-2466 1352

BRANCH OFFICES:

Bangalore: #105, Bhimajyothi Colony, West of Chord Road, Basaveshware Nagar, Above Coffee Day, Bangalore - 560 079 Tel.: +91-80-2323 0923 / 21
Fax: +91-80-2323 0924
Lucknow: 403, 4th Floor, Sahara Shopping Complex, Faizabad Road, Indira Nagar, Lucknow - 226 016. Tel.: +91-522-2346 535 / 2346 496 Fax: +91-522-23488 847
Hyderabad: 6-3-1090/1/, Uma Hyderabad House, 2nd Floor, Somajiguda, Raj Bhavan Road, Hyderabad - 500 082. Tel.: +91-40-6550 4133 / 2331 3300
Fax: +91-40-2331 3312
Gurgaon: Plot No. 372, Uday Vihar Phase II, Gurgaon - 122 015, Hararyana.
Tel.: +91-124-2843 300 / 2843 600 Fax: +91-124-2843 399
Tiruvanthapuram: Megha Complex, Ground Floor, TC No. 26/857, Women College Road, Thycaud, Thruvanthapuram - 695 014
Tel.: +91-471-2324 064 / 2320 082 Fax: +91-471-2320 082
Vadodara: 401/402, Manan Commercial Landmark, Opp. Dhara Hospital, Above HDFC Bank, Gotri Road, Vadodara - 390 015 Tel.: +91-265-2342 929
Fax: +91-265-2352 924
Pune: Kailaschandra 128 / 1A, 2nd Floor, Paud Road, Kothrud, Pune - 411 038
Tel.: 098220 43622
Chandigarh: S.C.O. 208-209, 4th Floor, Basera Building, Sector-34 A, Chandigarh - 160 022, U.T. Tel.: +91-172-4090 001-07 / 4090 009
Fax: +91-172-4090 008



Striving to become the best individuals, we endeavour to foster the best team. Performing sensibly, we try to achieve the best efficiency. Working innovatively, we seek to make the best products. Listening patiently, we excel to offer the best service. So, no matter what your needs are, come to us, GET THE BEST

LABINDIA reserve the right to change specification without notice as part of its continuous programme of product development.